

Oct 07, 2017 Version 1

, , ,

Ischemia-reperfusion-induced acute kidney injury V.1

PLOS One

DOI

dx.doi.org/10.17504/protocols.io.j8gcrtw

Laura Ruiz Remolina<sup>1</sup>

<sup>1</sup>University of Salamanca (USAL) Spain



Laura Ruiz Remolina

## OPEN ACCESS



DOI: dx.doi.org/10.17504/protocols.io.j8gcrtw

External link: https://doi.org/10.1371/journal.pone.0188204

Protocol Citation: Laura Ruiz Remolina 2017. Ischemia-reperfusion-induced acute kidney injury. protocols.io

https://dx.doi.org/10.17504/protocols.io.j8gcrtw

## Manuscript citation:

Ruiz-Remolina L, Ollauri-Ibáñez C, Pérez-Roque L, Núñez-Gómez E, Pérez-Barriocanal F, López-Novoa JM, Pericacho M, Rodríguez-Barbero A (2017) Circulating soluble endoglin modifies the inflammatory response in mice. PLoS ONE 12(11): e0188204. doi: 10.1371/journal.pone.0188204

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: October 07, 2017

Last Modified: January 15, 2018

Protocol Integer ID: 8168

Keywords: Ischemia-reperfusion, acute kidney injury



## Materials

## **MATERIALS**

- X NaCl 0.9%
- **⋈** Buprenorphin **Catalog #**961425
- Surgical Staples Catalog #138007
  ■
- xurgical material
- **23**G Needles **Catalog #**4657667
- **S** Gauzes
- X Artery Clamp



Once the mouse is anesthetized, we place it face up on the electric blanket, to maintain the body heat.

1

We make with scissors an incision in linea alba, at the level of the kidneys. The incision will have to be as small as possible, but of sufficient size to work comfortably.

2

Trying to manipulate the intestinal package as little as possible, we locate the renal pedicel (located behind the intestines, a deep red color). We will do the ischemia in the left kidney of the mouse (during the operation, the one on our right).

3

Clean the conjunctive around the pedicle with a pair of forceps and clamp it with the artery clamps, thus performing ischemia. to. a. If the ischemia is well made the kidney will darken, becoming black in a few minutes. b. In older mice the technique is complicated because of the greater amount of fat. c. Once the ischemia was performed, the analgesic BUPREX was injected with the mouse: 200-300ul dilution 1:20 (in saline). Ex: 15ul BUPREX + 300ul NaCl.

4

We maintain ischemia for 30 minutes, keeping the mouse covered with gauze and watching at all times for no bleeding.

5



When removing the artery clamps that are producing ischemia, we see how it reperfuses the kidney in just a few seconds.

6

The peritoneum is continuously sewed, azole powders are added and the skin is closed with staples.

7

We injected 200ul of saline to the mouse so that it is hydrated.

8