


Apr 11, 2022 Version 2

Western blotting in *Chlamydomonas reinhardtii* V.2

 PLOS One

DOI

dx.doi.org/10.17504/protocols.io.14egnjnpv5dy/v2

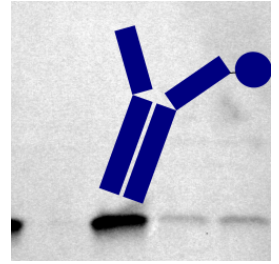

João Vitor Molino¹

¹University of São Paulo



Joao Vitor Molino

Ronin Institute, University of California, San Diego

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.14egnjnpv5dy/v2

External link: <https://doi.org/10.1371/journal.pone.0192433>

Protocol Citation: João Vitor Molino 2022. Western blotting in *Chlamydomonas reinhardtii*. [protocols.io](https://dx.doi.org/10.17504/protocols.io.14egnjnpv5dy/v2) <https://dx.doi.org/10.17504/protocols.io.14egnjnpv5dy/v2> Version created by [Joao Vitor Molino](#)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 11, 2022

Last Modified: April 11, 2022

Protocol Integer ID: 60613

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

This protocols describe the steps to perform a western blot in *Chlamydomonas reinhardtii* cell lysate and supernatant samples.

Guidelines

Prepare cell material by cultivating in liquid media.

Before start

- Check antibody dilutions
- Check buffers disponibility
- Prepare 5% milk solution



Sample preparation - Supernatant

1. Centrifuge algae culture at 2000xg for 10 min
2. Recover supernatant

00:10:00 Centrifugation time

Sample preparation - Lysate

- 2 1. Centrifuge algae culture at 2000xg for 10 min
2. Remove supernatant, and re-suspend cells in lysis buffer (50 mM Tris·HCL (pH 8.0), 0.1% Triton X-100), concentrating cells 100-fold.

00:10:00 Centrifugation time

Sonication

- 3 1. Sonicate using appropriate sonication tip. Duty cycle: 0.5 | Cycle: 1.0 s | Amplitude: 20% | Duration: 30 s
2. Centrifuge for 15 min at 20000xg to remove cells debris and recover soluble proteins
3. Quantificate soluble protein

00:15:00 Centrifugation time

Gel electrophoresis | SDS-PAGE

- 4 1. Load 30 µg of total soluble protein (TSP) per lane in a 12% SDS-PAGE
2. Transfer proteins to a nitrocellulose membrane
3. Block the membrane with 5% milk solution
4. Probe the desired protein with the specific antibody, diluted in 5% milk solution
5. Wash 2 times with TBST (0.2 M Tris, 1.37 M NaCl, 0.1% Tween-20, pH 7.6)
6. Add secondary antibody if required

30 µg total soluble protein per lane